

Streptavidin, HRP-Labeled

molecular biology grade



Catalog No.
474-3000

Size
1 mL

DESCRIPTION

Streptavidin is a 60,000 dalton protein isolated from the bacterium *Streptomyces avidinii*. Streptavidin has been shown to bind four molecules of biotin with high affinity ($K_D=10^{-14}$ M) (1). Electrophoretically pure streptavidin is linked to horseradish peroxidase using a modification of the periodate method of Nakane and Kawaoi (2).

FORM

The conjugate is provided in a liquid buffer containing a proprietary stabilizer and an anti-bacterial agent. It is prepared with molecular biology grade chemicals and dispensed into RNase/DNase-free sterile vials.

STORAGE/STABILITY

Store at 2-8°C. Stable for a minimum of 1 year from date of receipt at 2-8°C as an undiluted liquid. Dilute immediately before use.

ENZYME:PROTEIN RATIO

Molar peroxidase:streptavidin ratio = 2.5:1.

CONCENTRATION

The concentration of streptavidin is 0.1 mg/mL in a 1.0 mL volume.

APPLICATIONS

Peroxidase labeled streptavidin is suitable for use in Northern blotting, Southern blotting, plaque and colony hybridizations, *in situ* hybridization, and immunohistochemistry (brief protocols described below) applications. This conjugate may also be used for ELISA and immunoblotting procedures (See References 3-6).

SUGGESTED WORKING DILUTIONS

Different assay conditions require that serial dilutions of all reagents be performed to determine optimal working concentrations. Prepare the working dilution immediately before use. Long-term storage at a working dilution may result in enzyme inactivation and performance loss. Do not use sodium azide in the diluent. For suggested starting dilutions, see the appropriate protocol.

SUGGESTED PROTOCOLS

All steps are at room temperature unless otherwise noted.

Southern Blotting, Northern Blotting, Plaque and Colony Hybridizations

Following hybridization with a biotinylated probe and post-hybridization washing:

1. Place membrane in a small container or hybridization bag, block with KPL's Detector Block (See RELATED PRODUCTS), or other appropriate blocking solution, for 30 minutes.
2. Dilute peroxidase-labeled streptavidin 1:100-1:1000 in blocking solution, using at least 0.25 mL per cm² membrane. Incubate membrane for 20 minutes.
3. Transfer membrane to a clean container and wash with KPL's Biotin Wash Solution (See RELATED PRODUCTS). Wash 3 times for 5 minutes each using at least 0.4 mL wash solution per cm² membrane.
4. Detect using TMB Membrane Substrate, LumiGLO® Chemiluminescent Substrate (See RELATED PRODUCTS) or other appropriate peroxidase substrate following manufacturer's protocols.

In Situ Hybridization

Following hybridization of tissue or cells with a biotinylated probe and post-hybridization washing:

1. Dilute peroxidase-labeled streptavidin 1:20-1:200 in an appropriate diluent and apply approximately 100 µL to specimen. Cover slide to prevent evaporation. Incubate in a 37°C humidified chamber for 20 minutes. The optimal dilution of HRP-SA must be determined experimentally.
2. Immerse slide in a coplin jar containing KPL's Biotin Wash Solution (See RELATED PRODUCTS), or other appropriate wash solution. Wash 3 times for 5 minutes each.
3. Detect using TrueBlue® (See RELATED PRODUCTS) or other appropriate peroxidase substrate following manufacturer's instructions.
4. Apply a red counterstain, such as Orcein (See RELATED PRODUCTS), according to manufacturer's instructions.
5. Dehydrate and mount the slide with an organic mounting media such as Permount.

Immunohistochemistry

Following incubation of the specimen with primary and biotin-labeled secondary antibody:

1. Dilute peroxidase-labeled streptavidin 1:10-1:100 in 10% Normal Goat Serum (See RELATED PRODUCTS). Flood the slide with diluted peroxidase-labeled streptavidin. Incubate for 30 minutes.
2. Soak the slide in PBS or Tris-HCl for 5 minutes.

3. Detect using TrueBlue® (See RELATED PRODUCTS) or other appropriate peroxidase substrate, and counterstain if desired.

PRODUCT SAFETY AND HANDLING

This product is considered non-hazardous as defined by The Hazard Communication Standard (29 CFR 1910.1200). Avoid contact with skin and eyes. In case of contact or spillage, clean with copious amounts of water. Product may be disposed via a sanitary sewer.

REFERENCES

1. Holmberg, A., et al. (2005). *Electrophoresis*, Feb;26(3):501-10.
2. Nakane, P.K. and Kawaoi, A. (1974). *J. Histochem. Cytochem.* 22: 1084-1091.
3. Brigati, D.J., et al. (1983) *Virology*, 126: 32-50.
4. Harlow, E. and Lane, D. (1988) Antibodies: A Laboratory Manual. Cold Spring Harbor Laboratory Press, NY.
5. Sambrook, J., et al. (1989) Molecular Cloning: A Laboratory Manual. Cold Spring Harbor Laboratory Press, NY.
6. Crowther, J.R. (1995) Methods in Molecular Biology: ELISA: Theory and Practice. Vol. 42, Humana Press, NJ.

RELATED PRODUCTS

Detector Block	Cat. No. 71-83-00
Biotin Wash Solution (10X)	Cat. No. 50-63-06
10% Normal Goat Serum	Cat. No. 71-00-27
LumiGLO® Chemiluminescent Substrate	Cat. No. 54-61-02
TMB Membrane Substrate	Cat. No. 50-77-18
Orcein Counterstain	Cat. No. 71-01-01
TrueBlue®	Cat. No. 50-78-02

Visit www.kpl.com for a complete list of KPL Products.

The product listed herein is for research use only and is not intended for use in human or clinical diagnosis. Nothing disclosed herein is to be construed as a recommendation to use this product in violation of any patents. The information presented above is believed to be accurate. However, said information and product are offered without warranty or guarantee since the ultimate conditions of use and the variability of the materials treated are beyond our control. We cannot be responsible for patent infringements or other violations that may occur with the use of this product. No claims beyond replacement of unacceptable material or refund of purchase price shall be allowed. All claims regarding product performance must be made within 30 days following date of delivery.

Limited Use License

The purchase of this product conveys to the buyer the non-transferable right to use the product in research conducted by the buyer. The buyer cannot sell or otherwise transfer this product or materials made by use of this product to a third party or otherwise use this product or materials made with this product for Commercial Purposes without written approval of KPL, Inc. For additional information, please visit www.kpl.com or the KPL catalog. To obtain a license or approval to use this product for purposes other than those permitted above, contact Senior Director of Sales and Marketing at (301) 948-7755.