

Protein G Agarose Kit

Catalog No.
553-51-00

**Kirkegaard & Perry
Laboratories**

KPL

INTRODUCTION

Protein G Agarose consists of recombinant protein G, which is produced in *E. coli* and, after purification, is covalently immobilized onto 4% cross-linked agarose beads. Protein G agarose is suitable for the isolation of IgG antibodies using column or immunoprecipitation methods. DNA sequencing of native protein G (from Streptococcal group G) has revealed two IgG-binding domains as well as sites for albumin and cell surface binding (1-6). Protein G has been designed to eliminate the albumin and cell surface binding domains to reduce nonspecific binding while maintaining efficient binding of the Fc region of IgGs. With the removal of these binding domains, Protein G can be used to separate albumin from crude human IgG samples (7).

Covalently coupled Protein G Agarose has been widely used for the isolation of a wide variety of immunoglobulin molecules from several mammalian species. Protein G has greater affinity for many more mammalian IgGs than Protein A (Table 1).

CONTENTS

The following components are found in the kit:

Protein G Agarose	Cat. No. 223-51-00
Binding/Wash Buffer, 5X Conc.	Cat. No. 50-70-01
Elution Buffer, 10X Conc.	Cat. No. 50-68-01
Storage Buffer	Cat. No. 50-69-01
Disposable Columns (2)	Cat. No. 80-00-10

- Protein G Agarose is supplied in a total volume of 15 ml consisting of 10 ml Protein G Agarose suspended in 20% ethanol/PBS.
- Binding/Wash Buffer is a 5X concentrate consisting of 0.5 M Sodium Phosphate and 0.75 M NaCl, pH 7.4.
- Elution Buffer is a 10X concentrate containing 2 M Glycine, pH 2.85.
- Storage Buffer is ready to use at 0.01 M NaH₂PO₄, 0.15M NaCl, 2.7 mM KCl, pH 7.4, 20% ethanol.

Also provided are 2 empty disposable columns with two sintered polyethylene frits with a pore size of 50-150 µm and reusable caps. The frits protect the agarose from running dry under gravitational buffer flow.

STORAGE/STABILITY

Store at 2-8°C. Stable for a minimum of 1 year from date of receipt when stored at 2-8°C. Non-sterile.

Note: Storage of the binding/wash buffer concentrate at 2-8°C may result in the appearance of salt crystals due to decreased solubility at reduced temperatures. Before preparing the 1X working solution, warm the binding/wash buffer at 37°C until all crystals have dissolved. Mix well by swirling vigorously, then proceed as described below. Once redissolved, this will have no effect on buffer performance.

Table 1. Relative Affinity of Immobilized Protein G and Protein A for Various Antibody Species and Subclasses of polyclonal and monoclonal IgGs (8).

Species/Subclass	Protein G	Protein A
MONOCLONAL		
Human		
IgG ₁	++++	++++
IgG ₂	++++	++++
IgG ₃	++++	---
IgG ₄	++++	++++
Mouse		
IgG ₁	++++	+
IgG _{2a}	++++	++++
IgG _{2b}	+++	+++
IgG ₃	+++	++
Rat		
IgG ₁	+	---
IgG _{2a}	++++	---
IgG _{2b}	++	---
IgG _{2c}	++	+
POLYCLONAL		
Rabbit	+++	++++
Cow	++++	++
Horse	++++	++
Goat	++	---
Guinea pig	++	++++
Sheep	++	+/-
Pig	+++	+++
Rat	++	+/-
Mouse	++	++
Chicken	+	---
Human IgG	++++	++++
Human IgM	+	---
Human IgE	+	---
Human IgA	+	---

--- (weak or no binding) → ++++ (Strong binding)

SPECIFICATIONS

Ligand density:

~2 mg Protein G/ml

	drained gel
Bead structure:	4% highly cross-linked agarose
Bead size range:	45 - 165 μ m
Recommended working pH:	3 - 9
Binding capacity:	\geq 18mg/ml Human IgG

Note: Different immunoglobulins derived from the same species and from the same subclass can demonstrate deviations in the binding capacity; Protein G may hydrolyze at low pH.

PROCEDURE

PURIFICATION OF IgG MOLECULES

1. **Buffer Preparation:** Prepare all buffers necessary for this procedure prior to starting.
 - a. **Wash/Binding Buffer:** Dilute the wash/binding buffer 1:5 in dH₂O in a clean container (e.g. 20 ml buffer concentrate + 80 ml dH₂O). Prepare enough diluted buffer based on the sample quantity to be dialyzed and ~30X the quantity of agarose used for purification. Store prepared buffer up to 1 week at 2-8°C. Final 1X concentration of buffer is 0.1 M Sodium phosphate, 0.15 M NaCl, pH 7.4.
 - b. **Elution Buffer:** Dilute the elution buffer 1:10 in dH₂O in a clean container (e.g. 10 ml buffer concentrate + 90 ml dH₂O). Prepare enough diluted buffer based on 4X the quantity of agarose used. Store prepared buffer up to 1 week at 2-8°C. Final 1X buffer concentration is 0.2 M Glycine, pH 2.85.
 - c. **Storage Buffer:** Ready to use.
2. **Sample Preparation:** To insure proper ionic strength and pH are maintained for optimal binding, it is necessary to dilute serum samples, ascites fluid or tissue culture supernatant at least 1:1 with binding buffer. Alternatively, the sample may be dialyzed overnight against wash/binding buffer. KPL recommends using a 12,000 MW cutoff dialysis tubing with at least two buffer exchanges. Remove any particulate matter from the sample by centrifugation or filtration through a 0.8 μ m filter.
3. **Column and Resin Preparation:**
 - a. Pour 20% Ethanol in the bottom of a petri dish or in a flat bottom container. Float the frit on top of the ethanol. Using the large round end of a 1 ml pipet tip, press the frit firmly into the ethanol to force air out. Repeat this step until the frit is completely wet.
 - b. Push the frit into the barrel of the column until it rests firmly on the bottom.
 - c. With the cap removed, clip the end of the column to create a hole to allow liquid to flow through.
 - d. Wash the frit with 5 column volumes of 1X Wash/Binding Buffer.

- e. Prepare a 1:1 suspension of resin in 1X Wash/Binding buffer. The required amount of agarose per mg immunoglobulin to be purified can be estimated by the binding capacity.

Recommended Column Volumes

Antibody Source	Recommended bed volume (ml) per ml sample
Immune Serum	2 ml
Tissue Culture Supernatant (with 10% fetal bovine serum)	0.2 ml
Tissue Culture Supernatant (serum-free)	0.01 ml
Ascites Fluid	2 ml

- f. Pour slurry into column. Allow column to flow by gravity to pack the column bed.
- g. Equilibrate the packed affinity resin with 10 column volumes (CV) of the wash/binding buffer. For example, if the packed bed is 1 ml, equilibrate with 10 ml Wash/Binding Buffer.
4. **Sample Purification**
 - a. Gently apply sample to the column by layering onto the top of the resin. Be careful not to disturb the bed surface.
 - b. Wash column with 10 CV of the 1X Wash/Binding buffer, or until the absorbance of eluate at 280 nm approaches the background level.
 - c. Before beginning the elution step, set up enough tubes to collect the entire elution volume as 1 ml fractions (4 CV will be used to elute the antibody). To each collection tube add 240 μ l 5X Wash/Binding Buffer. To elute the antibody, gently add 1 ml 1X Elution Buffer to the top of the resin collecting the eluate in a prepared collection tube. Repeat until the entire volume has been collected up to 4 column volumes. **Note:** If the eluate is to be collected in a single bulk volume, add 240 μ l 5X Wash/Binding Buffer per ml Elution Buffer to the collection vessel before starting the elution. Elution of bound immunoglobulin can be monitored by absorbance at 280 nm, if desired.
5. **Column Regeneration:** Once the sample has been eluted, wash the affinity matrix with 2 CV of elution buffer. Re-equilibrate the column with at least 10 CV of 1X Wash/Binding Buffer. When column is equilibrated, pH of eluate will be the same as that of the wash/binding buffer.
6. **Clean-in-Place:** With certain applications, substances which contain denatured proteins or lipids do not elute in the regeneration procedure. The following steps can be taken to clean the column:

- a. To remove strongly bound hydrophobic proteins, lipoproteins and lipids, wash the column with a non-ionic detergent (e.g. 0.1% Triton X-100) at 37°C, with a contact time of ~1 minute.
 - b. Immediately re-equilibrate the column with 5 - 10 CV of 1X Wash/Binding Buffer.
 - c. As an alternative, wash the column with 70% ethanol. Allow the column to stand for 12 hours. Re-equilibrate the column with 5 - 10 CV of 1X Wash/Binding Buffer.
 - d. To remove precipitated or denatured substances, wash the column with 2 CV of 6M guanidine hydrochloride. Immediately re-equilibrate the column with 5 - 10 CV of 1X Wash/Binding Buffer (see step 5).
6. **Resin Storage:** Store affinity matrix in Storage Buffer at 2-8°C. **Do not** store the matrix frozen or at room temperature. The matrix can be stored in the column by sealing the outlets or removed from the column and stored as a slurry.

PRODUCT SAFETY AND HANDLING

This product is considered non-hazardous as defined by The Hazard Communication Standard (29 CFR 1910.1200). Avoid contact with skin and eyes. Consult the Material Safety Data Sheet for precautions and disposal instructions.

3. Sjöbring, U., et. al. (1988). *J. Immunol.*, 140: 1595 - 1599.
4. Fahnestock, S. (1987). *TIBS*, 79 - 83.
5. Olsson, A., et. al. (1987). *Eur. J. Biochem.* 168: 391 - 324.
6. Fahnestock, S., et. al. (1986). *J. Bacteriol.*, 176: 870 - 880.
7. Goward, C.R., et. al. (1990). *Biochem. J.*, 267: 171-177.
8. Harlow, E. and Lane, D. eds.(1988). Antibodies, A Laboratory Manual. Cold Spring Harbor Laboratory, N.Y., 617-618.

RELATED PRODUCTS

Protein G Agarose	Cat. No. 223-51-01
Protein A Agarose Kit	Cat. No. 553-50-00
Protein A Agarose	Cat. No. 223-50-01

See KPL's catalog for a complete list of antibodies, western blotting and ELISA kits.

The product listed herein is for research use only and is not intended for use in human or clinical diagnosis. Nothing disclosed herein is to be construed as a recommendation to use this product in violation of any patents. However, said information and product are offered without warranty or guarantee since the ultimate conditions of use and the variability of the materials treated are beyond our control. We cannot be responsible for patent infringements or other violations that may occur with the use of this product. No claims beyond replacement of unacceptable material or refund of purchase price shall be allowed.

L-287-03

REFERENCES

1. Guss, B., et.al. (1986). *EMBO J.*, 5: 1567 - 1575.
2. Akerstrom, B., Nielsen, E. and Bjorck, L. (1987). *J. Biol. Chem.*, 262: 13388 - 13391.